

---

# OPERATION MANUAL

---

## *Model LET-2*

June 2004



**Far West Technology, Inc.**  
330 South Kellogg Ave, Suite D  
Goleta, CA 93117

## GENERAL INFORMATION

---

This instrument is manufactured in the United States of America by:

Far West Technology, Inc.  
330 South Kellogg Ave, Suite D  
Goleta, CA 93117-3814

Tel: 805.964.3615

Fax: 805.964.3162

Email: [support@fwt.com](mailto:support@fwt.com)

Web Site: [www.fwt.com](http://www.fwt.com)

Far West Technology has been manufacturing radiation measuring devices since 1972.

## REPAIR SERVICE

---

Although we design and manufacture our instruments to a high standard, we realize that repairs are sometimes necessary. If you believe service is needed on this instrument please call our service department before shipping the instrument to us for repair; often we can help you with simple problems. If you do decide to return it to us for repair then please include:

1. Contact person's name
2. Organization or Company name
3. Address
4. Phone number of contact person
5. Description of the problem
6. Anything else you may think important

We will inform you of the repair charges and wait for your authorization before we repair your instrument.

## TABLE OF CONTENTS

---

I. INTRODUCTION .....	2
II. PHYSICAL CHARACTERISTICS OF THE INSTRUMENT .....	2
III. FILLING THE INSTRUMENT WITH GAS .....	3
IV. ELECTRICAL CHARACTERISTICS .....	3
V. INTERNAL ALPHA SOURCE .....	4
VI. TYPICAL SPECTRUM FROM <sup>252</sup> CF NEUTRONS.....	4
VII. TYPICAL OPERATION OF THE INSTRUMENT .....	5
VIII. DATA REDUCTION METHODS .....	5
IX. REFERENCES .....	5

## I. INTRODUCTION

The instrument, as shown in Figure 1, is a spherical tissue equivalent proportional counter. It is usually filled with tissue equivalent gas at a reduced pressure. An aluminum shell is used as a vacuum tight container for the sphere and is mounted on an aluminum plate, which holds the electrical and gas connectors.

The instrument is generally used to accumulate a pulse height spectrum proportional to the energy deposited in the sensitive volume. This spectrum may then be transformed into a distribution of absorbed dose in LET with the aid of computer processing (or, at less convenience, by hand).

Simple electronic processing of the pulse height data as it is accumulated can also be used to allow measurement of exposures in Rads-tissue directly. This technique involves scaling of the oscillator pulses in the common Wilkinson type ADC used in most analyzers.

## II. PHYSICAL CHARACTERISTICS OF THE INSTRUMENT

The detector is a spherical cavity in tissue equivalent plastic (Shonka Type A-150)<sup>1,2</sup> with a 2.24 inch (5.69 cm) internal diameter. An aluminum can surrounds the TE plastic that provides electrostatic shielding and serves as a vacuum tight container. The aluminum is 0.063 inch (0.159 cm) thick. The plastic sphere is 0.125 inches (0.318cm) thick, which is sufficient for proton equilibrium at least to 20 MeV.

The collecting wire is positioned on a diameter of the sphere and is surrounded by a field equalizing helix. This design is similar to those originally published by H. H. Rossi and associates.<sup>6-8</sup> The collecting wire is .005 inches (0.0127cm) in diameter. The inside diameter of the helix is 0.125 inches (0.318cm).

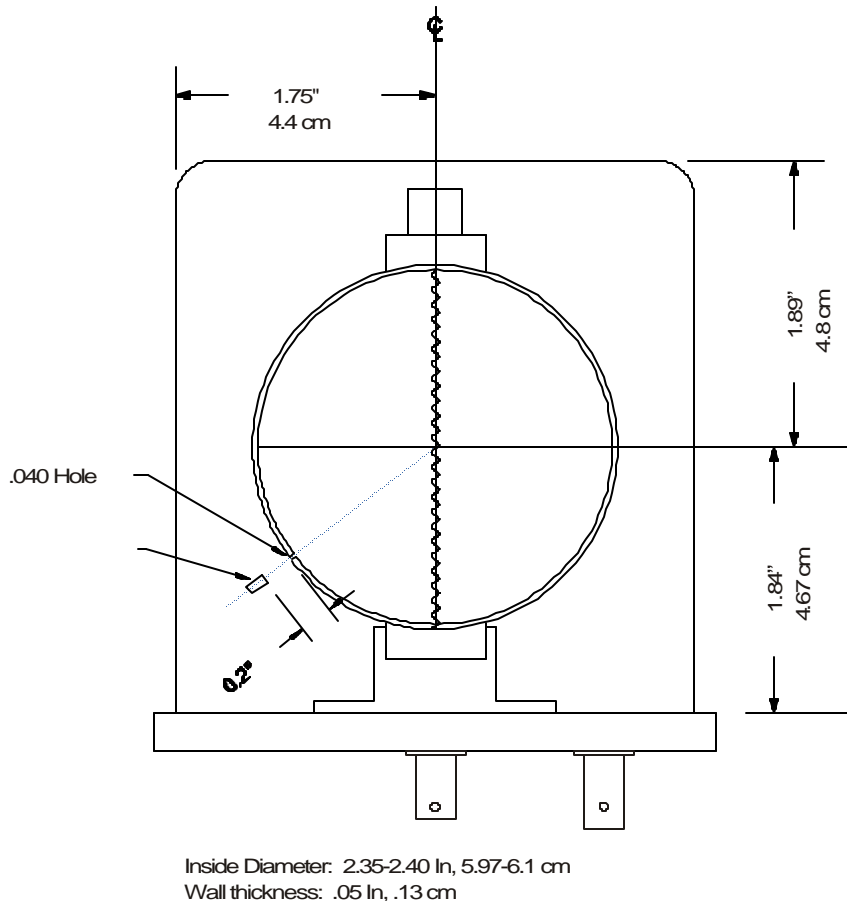


Figure 1 Dimensions of LET-2 Detector

### III. FILLING THE INSTRUMENT WITH GAS

---

The standard instrument is provided with a Quick-Connect manufactured by Crawford Fitting Company. This connector mates with a Swagelok B-QC4-D-400 DESO-type connector. These connectors contain spring loaded plungers that seat against elastomeric o-rings providing a gas-tight seal. Experience will allow the gas filling tube to be "snapped off" the instrument with no detectable change in instrument gain. Dirt must not be allowed to get into the Quick-Connect. It will invariably cause instrument gain shifts with time. When a bellows valve is used, this is not a problem.

The usual procedure for gas filling is, initially, a pump down to 5 to 10 microns with a good quality rotary pump. Liquid nitrogen cold traps may be useful, but have not been found necessary. The instrument is then filled to about 50 cm Hg with tissue equivalent gas and pumped down to 5 to 10 microns. This procedure may be repeated if the instrument has not been in use for some time.

The instrument is next filled to the proper pressure for operation. This is determined by the tissue equivalent gas density. Our usual gas mixture is 29.9% CO<sub>2</sub>, 2.75% N<sub>2</sub> and 67.54% CH<sub>4</sub> by volume with a density of 1.062 grams per liter at 20 c and 760 mm Hg. The instrument cavity is 2.24 inch in diameter (5.69 cm) and thus a pressure of 1.26 cm Hg will simulate a cavity of about  $1 \times 10^{-6}$  meter diameter in density 1.00 tissue. This effective diameter can be made smaller or larger by variation of the gas pressure. A more recent tissue equivalent gas mixture<sup>5</sup> composed of 39.6% CO<sub>2</sub>, 5.4% N<sub>2</sub> and 55% propane has been used with equivalent results. Addition of 10% isobutane will provide better operation at very high multiplications.

Gas purity is of considerable importance in counter operation. Poor resolution, gain or a low voltage arc-over point are evidences of gas problems. Water in the gas will invariably cause arc-over. A dew point of -40°C or lower is recommended.

### IV. ELECTRICAL CHARACTERISTICS

---

A resistor divider network is located inside the aluminum housing. It provides proper distribution of voltages between the central electrode, helix and outer shell. In this design the central electrode is operated at a high positive potential. This voltage is applied to the connector marked HV and is normally +600 volts at a T. E. gas pressure of 1.48 cm Hg (2.0 micron effective diameter). The voltages needed for propane based T.E. gas operation are given below:

Counter Diameter	Gas Pressure	Operating Volts	Arc-Over Voltage
<i>Methane Base TE</i>			
1 micron	1.26 cm Hg	700	900
2 micron	2.52 cm Hg	850	1000
<i>Propane Base TE</i>			
1 micron	0.74 cm Hg	550	620
2 micron	1.48 cm Hg	600	680

Other gases and gas mixtures differ in Operating Volts and Arc-Over Voltage, and are typically lower because of higher gas gains. Arc-overs should be avoided due to the inevitable insulator carbon tracking which occurs. This can seriously degrade the chamber performance. Large arc-overs can transfer enough charge to destroy the input FET on some preamplifiers. If a solid-state preamplifier is used, it is usually necessary to limit the rate of rise of the high voltage so that switching transients do not destroy the first amplifier stage semi-conductor. The high voltage connector mates with an MHV cable connector (UG-932 A/U or equal). RG-59U or equivalent is satisfactory for the high voltage connection.

The signal from the detector appears on the connector block marked SIG. This connector mates with BNC cable connector UG-260/U. Low noise cable such as Microdot (Malco) 250/3834-0000 should be used for best low noise

performance. The cable length should be as short as possible to reduce input capacity. The detector capacity to ground is approximately 9 picofarads. The pulses produced are negative going.

The SIG connector should be connected to a low noise preamplifier. The type usually specified for cooled Ge(Li) detectors is suitable (Tennelec TC-174 or 175, or Ortec 142A or 142PC). Gain should be approximately 100 to 200 mV/MeV(Ge). The detector output is AC coupled so the preamplifier may not need an input coupling capacitor. This may result in lower noise. Removal of any protection diodes across the input FET will also lower the noise, but the rate of rise of the high voltage must be limited when turning it on or off, in this case. See Figure 2.

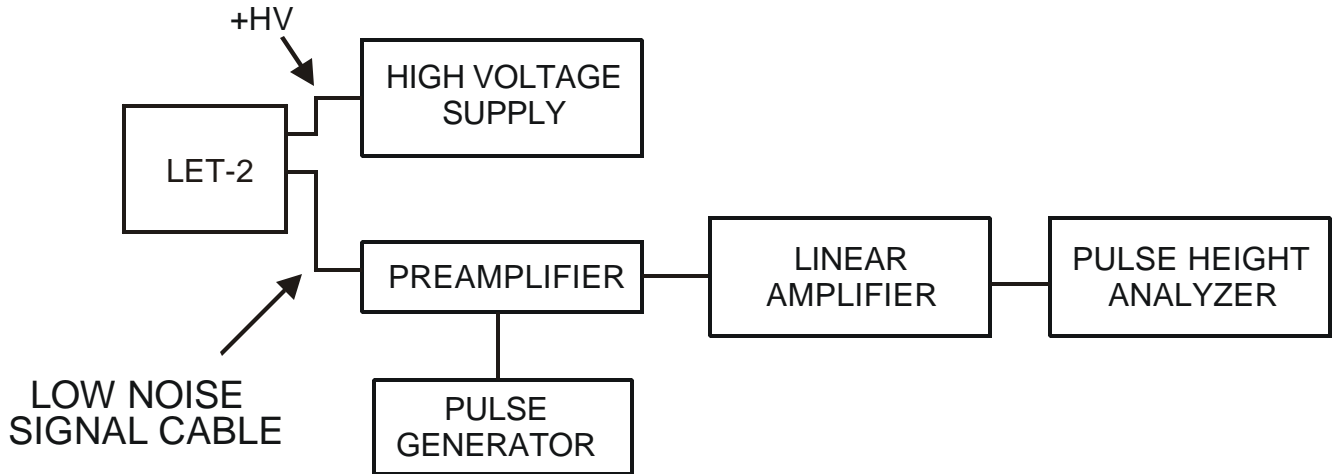


Figure 2 Electrical Interconnection Diagram for LET-2 Operation

## V. INTERNAL ALPHA SOURCE

The detector contains a  $^{244}\text{Cm}$  source for gain verification. The source is positioned so that the  $^{244}\text{Cm}$  alpha particles can enter the sphere through a collimator.  $^{244}\text{Cm}$  has a mean alpha energy of 5.80 MeV which averages to 81.72 keV/u over a range of 1u (1u =  $1 \times 10^{-6}$  meters) in tissue. Because of the collimator, the source produces a peak on a multichannel pulse height analyzer. The center of this distribution is characteristic of the LET of the alpha particle averaged over the detector diameter. Exact work can benefit from fitting a parabola to the upper half of this peak in order to find the precise position of the maximum. The alpha source resolution is relatively poor due to a compromise between source strength, collimator opening and useful count rate.

The standard LET-2 is supplied with a gravity operated source. It is activated by turning the detector on its side with the lettering on the baseplate upside down. If the LET-2 is supplied with the optional magnetically operated source, then there is a mark on the housing for the location of the supplied magnet. Place the magnet on the mark to activate the source.

A plot is included with each detector showing the alpha peak produced by its internal source. Data are included on the graph giving the high voltage used, gain settings of the linear amplifier, gain settings of the pulse height analyzer and the gas filling pressure.

## VI. TYPICAL SPECTRUM FROM $^{252}\text{CF}$ NEUTRONS

Plots of the pulse height spectrum derived from a standard  $^{252}\text{CF}$  exposure and from the internal  $^{244}\text{Cm}$  source are enclosed with each detector. These data were taken with standardized electronics as a quality control measure and are not representative of the best in low noise circuitry. The channel numbers written on the graph near the proton "edge" may be used to calculate resolution. These values so obtained are usually 10 to 11 percent.

## VII. TYPICAL OPERATION OF THE INSTRUMENT

---

The instrument is initially prepared as noted in the section "Filling of the Instrument with Gas", i.e., pump-down and flushing with T.E. gas, a second pump-down and filling to the required pressure. The Quick-Connect is snapped off freeing the instrument from the gas filling apparatus. The signal and high voltage cables are then connected (using low noise cable in the signal lead). The high voltage power supply is adjusted to provide +700 volts if the detector has an effective diameter of 1 micron. The signal from the preamplifier and through the linear amplifier is presented to a pulse height analyzer. A small magnet is placed on the counter where "magnet" is written to turn on the alpha source. The system gain is then adjusted to place the alpha peak in a convenient channel in the pulse height analyzer. For example, at an effective diameter of 1 micron, the average LET of the  $^{244}\text{Cm}$  alpha particles is 81.72 keV/u and if the peak is placed near channel 82, then the analyzer display is approximately 1 keV/u per channel.

After this preliminary work, an alpha spectrum is accumulated. Data should be accumulated for at least 10 minutes or more so that the position of the peak may be accurately determined. The alpha source is then turned off by removing the source magnet making the counter ready for data taking. Data are usually accumulated at several gain settings so that the limited dynamic range of the pulse height analyzer does not restrict the results. For example, the first run might be from about 40 to 400 keV/u (for the 400 channel analyzer). The second from 6 to 60 keV/u, the third from 1 to 10 and the fourth from noise up to 3 keV/u. This scheme allows overlap between the various segments so that they may be fitted together accurately. Obviously analyzer non-linearities must be known, as well as the precise gain shifts used to select the spectrum segments. We have used a sliding pulse generator to advantage in determining the analyzer non-linearities. A simple very stable mercury pulser can be used to measure the overall gain at the different settings.

After the data are accumulated it is good practice to take a noise spectrum for the highest gain segment. This will allow subtraction of the electronically produced noise and usually allows good data to be obtained down to about 0.5 keV/u. A second alpha spectrum should also be taken to establish detector drift characteristics, if any. Usually the detector can be expected to drift less than 1 channel out of 100 in 8 hours. Drift rates greater than this generally can be traced to gas leakage, either through pinholes or occasionally to the Quick-Connect. Since the latter is a o-ring sealed device a bit of dirt may allow some gas leakage. This can be cured in most cases by repeated operation of the Quick-Connect. The Quick-Connect leakage is not a problem if your detector is fitted with a bellows valve. Temperature changes can also cause gain shifts.

## VIII. DATA REDUCTION METHODS

---

Some methods of data reduction are discussed in References 6, 7, and 8. Invariably, computer aided processing is required because of the large number of data points gathered. Smoothing of the data before processing may also be helpful. Both linear and quadratic smoothing have been used with success.

If only certain segments of the data are needed initially a simple program can usually be written for one of the many programmable calculators now on the market. Such a program must include a smoothing routine if it is to be satisfactory.

## IX. REFERENCES

---

1. F. R. Shonka, J. E. Rose and G. Failla, "Conducting Plastic Equivalent to tissue, Air and Polystyrene", A/Conf. 15/p/753, Second United Nations International Conference on Peaceful Uses of Atomic Energy, June 1958.
2. J. B. Smathers, et al, October 1975, submitted to Medical Physics. A-150 composition given as H 10.2 +.1, C 76.8 +.5, O 5.9 +.2, N 3.6 +.2, Ca 1.8 +.1 and F 1.7 +.1 weight percent.
3. P. W. Benjamin, C. D. Kemshall and J. Redfern, "A High Resolution Spherical Proportional Counter", Nuclear Instruments and Methods, 59, 77 (1968).
4. B. Day, "A Portable LET Spectrometry System for the Calibration of the Concorde Radiation Meter", AWRE 057/69, Avg 1909

5. D. Srdoc, "Experimental Technique of Measurement of Microscopic Energy Distribution in Irradiated Matter Using Rossi Counters", *Radiation Research*, 43, 302 (1970).
6. H. H. Rossi and W. Rosenzweig, "A Device for the Measurement of Dose as a Function of Specific Ionization", *Radiology*, 64, 404 (1955).
7. H. H. Rossi and W. Rosenzweig, "Measurements of Neutron Dose as a Function of Linear Energy Transfer", *Radiation Research*, 2, 417 (1955).
8. H. H. Rossi, Chapter 2 in F. H. Attix and W. C. Roesch, *Radiation Dosimetry, Volume 1, Second Edition*, Academic Press, New York and London, 1968.
9. W. H. Grant, III, G. D. Oliver Jr, and B. A. Mitchel